VWR Bench top autoclaves

VAPOUR-Line *lite* II - VAPOUR-Line *lite* II P

INSTRUCTION MANUAL

European Catalogue Numbers:

481-0846 and 481-0847

Version: 3.0

Issued: 18th March 2019

Legal Address of Manufacturer

Europe

VWR International S.r.l.

Via San Giusto 85

20153 Milano (MI)

Italy

Tel. : 02-3320311/02-487791

<http://it.vwr.com>

Country of origin Italy

Table of Contents

Use of the manual	4
Safety information	4
Precautions during installation	4
Unit description	5
Electrical supply	6
Before using for the first time	6
Preliminary operations	7
Sterilisation temperature programming	7
Sterilisation time programming	7
Sterilisation cycle	8
Panel commands	8
Cleaning after use	8
General maintenance	9
Heating element replacement	11
Temperature/pressure rate	11
Spare parts	11
Setting the parameter safety key	12
Printer recorder (optional)	12
Safety precautions	13
Technical assistance	14
Warranty	14
Disposal	15

Use of the manual

Using this manual will help you use this instrument efficiently. The manual should be kept in a convenient place near the instrument for ease of use.

Safety information

- Use the device only for the indicated purposes
- The device must be correctly installed according to these instructions, before starting any operation
- Never disconnect the device by pulling the electrical cable from the socket
- Electrical cables should be replaced immediately when damaged
- While removing the item connected to a power source, pay attention to the electrical connection cable
- Always disconnect from the power source when:
 - A. Repairing or carrying out maintenance work; these operations must be carried out by qualified staff only
 - B. Cleaning the unit
 - C. Use original spare parts and accessories for any replacement

Do not use this device in the presence of explosive gases.

Precautions during installation

When the electrical system socket is not compatible with device, replace it paying attention to the following coded cable colours.

Brown: Live – Blue: Neutral – Yellow/green: Earth

Be aware that the voltage and frequency of the electrical system being used is compatible with the device.

The device must be connected to a ground wire. For any questions please contact a qualified electrician. Before starting up the device, be sure that the ventilation grids are totally free of obstruction.

Unit description

The steam steriliser 'VAPOUR-LINE *lite* II consists of:

Pressure gauge

The pressure gauge registers pressure (PSI). The pointer moves around the dial indicating precisely the pressure within the unit.

Pressure gauge adaptor

The pressure gauge adaptor connects the pressure gauge to the vessel cover.

Pressure regulator

The pressure regulator is a solid one-piece unit. Accurate pressure readings can be controlled and maintained by adjusting heat settings.

Air vent pipe

The pressure regulator fits over the vent pipe and allows excess pressure to be released.

Air vent cover lock

The air vent/cover locks automatically 'vents' or exhausts air from the vessel and acts as a visual indication of pressure in the unit.

Locking bracket

The locking bracket on the inside of the vessel body engages with the air vent/cover lock to prevent the cover from being opened when there is pressure in the unit.

Sealing ring

The sealing ring fits into the vessel cover and forms a pressure-tight seal between cover and body during sterilisation.

Overpressure plug

The overpressure plug is located in the vessel cover. It will automatically pop out and release steam in case the vent pipe becomes clogged and pressure cannot be released normally.

Litebox loading basket

Vessel and lid with two peripheral handles

Electronic device for thermoregulation

Main switch

Optional printer

Electrical supply

220 – 240 V – AC single phase, 1300 W.

This unit must be earthed. It is delivered with a mains plug with safety contacts.

If replacing the mains plug, the mains lead (coloured green and yellow) must be connected to the earth terminal on the new plug.

If using a mains supply without a protective earth, the unit must be connected to a safe earth.

Before using for the first time

- 1) Remove the air vent/cover lock from the vessel cover.

The cup portion of the air vent/cover lock, on the inside of the cover, unscrews from the pin.

- 2) Remove the sealing ring by simply pulling from the sealing ring groove.
- 3) Wash the cover and body with hot, soapy water. Do not submerge the cover in water or let the gauge come in contact with any liquid. The inside mechanism of the gauge may be damaged if it is submerged or if water is allowed to run over it. Scrub the sealing ring groove to remove the manufacturing oils. Scour the inside of the canner body and cover with a non alkaline, aluminium cleaner or a steel wool pad. Rinse with warm water and dry.
- 4) Re-insert the air vent/cover lock making sure the small gasket is in place.
- 5) Replace the sealing ring in the sealing ring groove.
The sealing ring is pre-lubricated; therefore, it is not necessary to apply oil to it. Frequent oiling may cause swelling of the sealing ring.
- 6) If the cover is difficult to open or close use a pastry brushes, a piece of cloth or your fingertips to apply a light coating of oil to the underside of the lugs on the vessel body. The cover should open and close easily.

Be sure the vent pipe is open before each use by holding the cover up to the light and looking through the vent pipe. If the vent pipe is blocked, excess pressure cannot be released through it.

Pressure will continue to build until the over pressure plug pops out of its cover opening.

Pressure is indicated on the dial gauge which registers the pressure inside the unit at all times. The pointer will move across the dial as pressure is increased or reduced within the vessel. The one-piece regulator is designed to rock only at 15 pounds pressure.

Do not strike the rim of the vessel body with any utensil as this will cause nicks which may damage the rim and allow steam to escape. Pouring water into a dry overheated vessel may crack the metal.

The vessel has been designed so the cover will go on in only one position and rotate only in the directions indicated for closing and opening. Press down on the cover handles to compress the sealing ring and make the cover easier to close. The cover has the words CLOSE/OPEN embossed on the top surface near a cover handle. Align the 'V' on the cover with the mark on the body handle

for proper cover and body alignment. Turn the cover in the direction indicated to close until the cover handles are centred directly above body handles.

Do not rotate the cover beyond this point.

When the cover is rotated to the closed position, the air vent/cover lock passes under the locking bracket. When the handles are aligned, one over the other, the pressure regulator is placed on the vent pipe and heat is applied, pressure will then begin to build within the vessel.

At this point, the air vent/cover lock lifts and locks the unit. The cover will remain locked as long as there is pressure in the vessel. When pressure is reduced completely, the air vent/cover lock drops allowing the unit to be opened.

Preliminary operations

- 1) This steam steriliser must only be used for a maximum pressure of 1.08 bar (121 °C).
- 2) The steam steriliser should be positioned in an area where there is adequate means for removal of steam.
- 3) Connect plug to mains supply.
- 4) Introduce distilled water or demineralised water into the sterilisation chamber to cover the heating system. Using distilled water avoids the production of calcareous deposits on the heating element. 20 minutes of sterilising will use approximately 500 ml of water.
- 5) Two displays are lit on time/temperature command panel.
- 6) Before programming, it is necessary to erase the protection of the parameters (safety keys), see paragraph 16.

Sterilisation temperature setting

- 1) Press 'SET' on lower display until Set.1 appears.
- 2) By using arrows increase or decrease the temperature. By keeping the button pressed, temperature will increase or decrease rapidly. For fine tuning, press the button once.
- 3) Press 'SET' to save to memory and repeat the same process for Set.2 parameters.

IMPORTANT. Set.1 AND Set.2 PARAMETERS SHOULD BE THE SAME VALUE

Sterilisation time setting

- 1) Press 'SET' until MA.t appears on the lower display..
- 2) Select the sterilization time using the arrows. The sterilisation time (in minutes) will be shown on the upper display.
- 3) Press 'SET' to save to memory.

Sterilisation cycle

- 1) Press the 'SET' button for 3 seconds to start the cycle. Actual temperature will be shown in the upper display (white), set temperature will be shown in the lower display (red)
- 2) When the sterilisation temperature has been reached, the time left to reach end of the cycle is shown on the lower display.
- 3) The end of the sterilisation cycle is indicated by an acoustic alarm for few seconds.
- 4) The cycle can be stopped at any time by simply pressing SET button for five seconds
- 5) At the end of vessel cooling, open the lid to take out the sterilised material.

Panel commands

SET	To start/stop the programme and change parameters to be set.
▼▲	To increase/decrease time and temperature
FNC	To enter service menu. A password will be required. To be used by service personnel.

NOTES:

All programming operations must be performed in quick sequence otherwise the programming device zeros. The programme can be stopped by pressing the 'SET' button for more than five seconds. The only parameters that need to be selected are (SP1) temperature and (t in 1) time. The selected parameters remain memorised even after the instrument is turned off.

Cleaning after use

- 1) The outside surface may be kept bright and shiny by cleaning with a good silver polish or simply by washing with soap and water. Iron and various minerals in water may darken the inside but this discoloration will in no way affect sterilisation. These stains are harmless and may be removed by using a solution of water and cream of tartar. For each quart of water, use one tablespoon of cream of tartar. Pour enough solution into the vessel to cover the discoloration (do not fill over 2/3 full), then close cover securely. Place regulator on vent pipe and heat until 15 pounds pressure is reached. Switch off and allow the unit to stand for two to three hours. Remove regulator, open and empty contents. Scour thoroughly with a soap impregnated steel wool cleaning pad; wash, rinse and dry.
- 2) Each time the vessel is washed, remove the sealing ring and wash in warm, soapy water, rinse, dry and replace in cover.

- 3) The air vent/cover lock may be removed for occasional cleaning or for replacing the small gasket. Wash all parts in warm, soapy water. Use a soft cloth or small nylon brush to clean the cover hole. The small gasket must be in place when re-assembling the air vent/cover lock.
- 4) The overpressure plug can be removed for cleaning by pulling it out of its opening from the bottom part of the cover. After cleaning, re-insert the plug by pushing the round side of the plug into the opening from the bottom part of the cover. When the overpressure plug is properly installed the word (TOP) will be visible when viewing the outside of the cover. The dented portion of the overpressure plug is visible when the bottom part of the cover is viewed. If the overpressure plug is ever forced out of its cover opening due to excess pressure while sterilising, it is important to have the vessel properly inspected before further use. Do not attempt to replace the overpressure plug and ask for servicing.
- 5) Be sure the vent pipe is open before each use. Clean the vent pipe with a small brush or pipe cleaner.
- 6) When your unit is not in use, invert the cover on the body and store in a dry place. Storing the unit with the cover locked on may deform the sealing ring.

General maintenance

- 1) CAUTION: To ensure safe operation and satisfactory performance, replace the overpressure plug every time you replace the sealing ring, or sooner if it becomes hard or deformed. In any case replace the overpressure plug every 2 years. Failure to follow these instructions could result in bodily injury or property damage.
- 2) If the body or cover handles become loose, tighten with a screwdriver.
- 3) If leakage of moisture or steam develops while using your unit, check the following possible causes:
 - The formation of a small amount of moisture under the pressure regulator is normal when heating first begins. This condensation is a result of the temperature of the pressure regulator being lower than the rest of the vessel. If excess condensation continues, the vent pipe may be loose and should be tightened with an adjustable wrench.
 - Leakage between the cover and body is usually caused by shrinkage of the sealing ring after prolonged use. Replace the sealing ring, overpressure plug and pressure gauge adaptor.
 - A small amount of leakage around the air vent/cover lock is normal when sterilising first begins. If leakage continues, the cover handles may not be fully aligned with the body handles and, therefore, the cover lock cannot engage. Clean the air vent/cover lock occasionally to ensure that it operates correctly.
 - A small amount of steam or moisture may be visible around the overpressure plug as sterilisation begins. This will stop when the overpressure plug seals. If leakage continues, clean or replace the overpressure plug.

- 4) If the pressure vessel becomes difficult to open or close and the sealing ring has expanded, replace the sealing ring.
- 5) **IMPORTANT:** In time, the sealing ring may shrink and the air vent/cover lock, small gasket, overpressure plug and pressure gauge adaptor may become hard and deformed. When this happens, replace the sealing ring, small gasket, overpressure plug and pressure gauge adaptor. The sealing ring, overpressure plug and pressure gauge adaptor should be replaced at least every 2 years.

Before inserting a new sealing ring, clean the sealing ring groove with a brush. To ensure safe operation of your unit, it is recommended that the pressure gauge adaptor and overpressure plug be replaced whenever the sealing ring is replaced.

- 6) **Anti-corrosion technique:** When two different metals, with different chemical properties, come in contact through an electrolyte an electron exchange is established. This process is known as galvanic corrosion which starts oxidation of the weakest metal. In an aluminium autoclave this natural effect can eventually lead to severe damage to the vessel.

To prevent this corrosion, all aluminium autoclaves are equipped with a special device called sacrificial body (Cat. No. 481-0721). This device is a small disc fixed with a screw to the vessel body and is made of a metal that is more easily oxidised than aluminium. This means that the galvanic current will flow through the sacrificial body which will be corroded instead of the autoclave vessel. If the sacrificial body has been completely worn out, the autoclave vessel will begin to corrode, so it is mandatory to periodically check the sacrificial body status and replace it before it wears out. It is advisable to order some spare sacrificial bodies in order to always run the autoclave with a sacrificial body in it.

Heating element replacement

Disconnect the electric supply, unscrew the two fixing nuts on the outside bottom of the sterilisation chamber and insert a new heating element. To re-assemble, repeat the same steps in reverse order. Ensure that the water level in the sterilisation chamber always covers the heating element.

Temperature/pressure rates

Temperature (°C)	100	102	105	109	112	115	119	120
Pressure (PSI)	0	1.45	2.90	5.80	7.25	10.15	13.05	14.50

Spare parts

	Description	Cat. No.
1	Manometer	481-0717
2	Heating element 200 V, 1300 W (110 V on request)	481-0777
3	Sealing ring	481-0719
4	Overpressure plug	481-0720
5	Main switch	481-0779
6	Sacrificial body	481-0721
7	Pressure regulator	481-0722
8	Controller for autoclave	481-0848
9	Printer	481-0840
10	Black start button	481-0839
	Printer paper	481-0775

Setting the parameter safety key

Selected parameters should be protected to avoid tampering:

- Press 'ENTER' several times until the word 'nnn' shows on the display
- Select a number other than 3 using the arrow keys and press 'ENTER'

This number will be requested any time the set parameter is changed.

The parameter protection is erased by doing the following:

- Press the 'ENTER' button several times until the word 'nnn' shows on the display
- Select number 3 using the arrows and press 'ENTER'

The programmer will now no longer request a password.

Printer recorder (optional)

The autoclave is supplied with a temperature printer connection.

Printer records:

- a. Date
- b. Batch number
- c. Temperature recorded during the segment of time programmed
- d. Sterilisation progressive number carried out automatically

How to insert the paper

Pressing the two levers (E), open the panel and insert paper as shown.

If the paper is correctly installed the 'pos F' LED light is on.

A flashing LED indicates that the paper is not properly installed.

By pressing button G you can scroll the paper.

Date:	2004/01/01	A
Name:	-----	
Batch:	1	B
96	10	
99	10	
105	10	
110	10	
115	10	
117	10	D
119	10	

Safety precautions

Opening and closing

The autoclave has a safety block when the lid is open. This block will prevent the lid falling on the hands of the operator. ALWAYS make sure that the block is working properly and that there's sufficient space between the instrument and the wall behind, in order to allow the lid to open completely and block to be activated.

After a cycle, the instrument should only be opened when temperature is below 70 °C and the inner pressure is at 0. Autoclaves have safety devices in order to avoid accidental opening. In this case DO NOT FORCE the instrument open.

Opening the autoclave when power is off is NOT recommended.

Personal Safety Systems (PSS)

All the safety devices on this autoclave will prevent accidental opening. However, sterilised loads can undergo thermal shock during a cycle that can cause breakages during opening, although in very rare circumstances.

The following PSS workwear should be near the instrument and always worn by personnel:

- Heat resistant gloves
- Heat resistant face mask (e.g. made of polycarbonate)
- Protective glasses
- Labcoat

Sterilisation of liquids

Please note that liquids in a load have a thermal capacity bigger than the steam in the chamber. For this reason they're always a "step behind" the temperature measured in the chamber. The more viscous or bulky the liquid is, the more this phenomenon may occur.

So, when the chamber is less than 100 °C, liquid inside the load can be at a temperature higher than this. Consequently if the autoclave is opened, liquid inside (that is at a much higher temperature) may come out in a turbulent manner and dirty the internal chamber or, (worst case) splash the operator. If the container is closed it can explode.

If you're sterilising liquids, always consider **at least** a 20 °C difference between chamber and temperature and load temperature. The instrument should therefore be opened at 50 °C.

We strongly recommend not to sterilise liquids in containers bigger than 2 litres.

Technical service

Web Resources

Visit the VWR's website at vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit.

vwr.com.

Warranty

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle end by handing it over to an authorised facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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VWR International AB
VWR International AG
VWR International Ltd