## Funktionsgenerator 250350

English  
manual  
at p. 7

2016.03.29 / HS

AC 250350

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2016.03.29 / HS

AC 250350

- 0.001 Hz to 10.0 MHz in one range
- Frequency control with smart, smoothly accelerating characteristic
- Step function: Jump through an overtone series easily – or through octaves – or decades (p. 9)
- Waveform: sine, triangle, square (bipolar) – square pulse, triangle pulse, ramp up & down (positive)
- Custom-defined waveform
- Sweep function – spot any resonances in a frequency range (p. 9)
- Built-in 10 W power amplifier

### QUICK GUIDE

**Preparation:** Turn down the *Output* knob almost completely. Turn the generator *On*.

**Waveform:** Select by pressing *Waveform*. Sine ( $\sim$ ) is good starting point.

**Frequency setting:** The *Frequency* knob is speed sensitive. Turn quickly to get near the desired frequency and slower to fine-tune. To go far, buttons  $\blacktriangle$  *Up* and  $\blacktriangledown$  *Down* may be used.

**Connections:** A speaker or vibrator normally connects to the *Power* jacks. Other devices connect to the *50 Ohm* output with coaxial cable. (The power output is disabled for frequencies exceeding 100 kHz.)

**Amplitude:** Adjust the level appropriately with the *Output* knob. The peak voltage for the output is indicated briefly in the display. The amplitude is also shown when the *Display* button is pressed.

**Step:** The fundamental frequency for the step function is set to the current frequency by pressing *f<sub>0</sub>/Mode* once.

Fast, repeated presses will switch between the modes harmonics (overtones), octaves and decades – indicated as *Harm*, *Oct* and *Dec*.

Use buttons  $\blacktriangle$  *Up* and  $\blacktriangledown$  *Down* to step through a series of overtones, octaves or decades.

**Sweep:** The first four times the button *Sweep* is pressed, a sweep parameter can be set. By the fifth press, the sweep starts. Yet another press stops the sweep and the generator is again in normal mode.

## Waveforms

The first three waveforms: sine, triangle and square – are the familiar signals with equal positive and negative extent.

The following four waveforms: square pulse, triangle pulse, ramp up and ramp down – has no negative values.

The display can show the signal amplitude, either when it is changed, or by pressing the *Display* button. In any case, the unloaded signal's amplitude – i.e. the positive peak value – is shown. Other popular specification of amplitude, such as effective value (RMS) or peak-to-peak (p-p), you will have to calculate yourself.

The last wave form: *Prog.* – is user defined. It can only be selected with the *Waveform* button if it is defined. (See section *Advanced use*).

## Frequency control

The *Frequency* knob is speed sensitive. A steady movement will give a frequency variation that is useful e.g. when searching for resonance frequencies in a column of air. Turning slower, 1 “click” on the knob will ultimately correspond to a change of the last digit by 1. And conversely, higher speeds result in greater and greater frequency changes.

We have chosen the acceleration characteristic of the knob after many practical experiments with common physics equipment. Such a choice will always reflect subjective preferences. Therefore it is possible to select other characteristics. This is described in the section *Advanced use*.

For large frequency changes, it can often be a shortcut to use the step function (see below).

The function generator 250350 is based on digital frequency synthesis, and therefore has a sharp upper cut-off frequency. Moreover, the analogue amplifier circuits impose their own limitations on the bandwidth. Sine waves are reproduced in the whole frequency range from 0.001 Hz to 10 MHz with very low distortion, while the other waveforms will have slightly rounded corners at high frequencies. This is a natural consequence of their harmonics being attenuated. However – even at 10 MHz, the diverse built-in waveforms remain recognizable.

When you define a waveform yourself (described in the section *Advanced use*), remember these limitations. You cannot make a 100 MHz function generator by defining a wave-form consisting of a sine wave that oscillates 10 times.

## Outputs

The *Power* output is used for power-consuming devices. This output operates in the frequency range from 0.001 Hz to 100 kHz. At higher frequencies the output is disabled.

Please note that the output is disabled completely in sweep mode even if only part of the sweep is above the limit of 100 kHz.

The *Power* output is thermally protected against overload, but should not be deliberately exposed to larger loads than specified (maximum output current: 1 A).

The *50 Ohm* output is active throughout the generator frequency range.

This output impedance is matched to 50  $\Omega$  coaxial cable, which is significant at high frequencies: If the cable is properly terminated with 50  $\Omega$ , there should be no reflections in the cable. In this setup, the amplitude of course decreases by half. Be aware that the display shows the amplitude “on the inside” of the output impedance, and therefore does not reflect the amplitude of the loaded output.

Note – check if the termination resistor can handle the deposited power (up to 0.5 W).

Normal use in the low frequency range (e.g. audio) does not require termination.

The *50 Ohm* output is short circuit protected, but you should not operate the unit for longer at full amplitude and high frequencies in a low impedance load, as this leads to substantial power dissipation within the generator.

The two previously mentioned outputs will have identical (adjustable) signal amplitudes when not loaded.

The *Sync Out* output is an extra output with fixed amplitude. The waveform of this output is always (positive) square pulses. Voltage levels are 0 and 5 V.

*Sync Out* is primarily intended as an external trigger signal to an oscilloscope. For small or noisy signals, it may prove helpful in order to get a stable image.

## The step function

In many applications you will study a physical system at a fundamental frequency and at different multiples of this frequency. The step function is created to facilitate this work.

The fundamental frequency for the step function is set to the current frequency by pressing  $f_0/Mode$  once. Fast, repeated presses will switch between the modes harmonics (overtones), octaves and decades – indicated as *Harm*, *Oct* and *Dec*.

The fundamental frequency is indicated in the display ( $f_0 = xxxx$ ).

The ▲ *Up* and ▼ *Down* buttons can then be used to jump up and down in the selected set of frequencies as described below. The current overtone, octave or decade is indicated by the display ( $N = xx$ ).

### Harmonic (Harm):

The frequency steps through the overtone series:  $f = f_0 \cdot N$ . Negative  $N$ -values should be interpreted as fractions: -2, -3 ... means  $1/2$ ,  $1/3$  ... ( $N$ -values 0 and -1 does not occur.)

### Octaves (Oct):

The frequency steps through octaves:  $f = f_0 \cdot 2^N$

### Decades (Dec):

The frequency steps through decades:  $f = f_0 \cdot 10^N$

## Application example – Step

In a study of standing waves on a string one finds a resonance at 175.0 Hz. The standing wave shows 5 anti-nodes (the string is 5 half wave-lengths long). Push  $f_0/Mode$  to set  $f_0 = 175.0$  Hz. If the display doesn't already display Harm, press again until it does. The display now shows  $N = 1$ .

Press ▼ *Down* four times to set  $N$  to -5 – the frequency is now  $f_0/5 = 35.00$  Hz.

This is the fundamental of the string and you observe a standing wave with 1 anti-node. In order to set the generator's fundamental frequency to 35.00 Hz also, press the  $f_0/Mode$  once. Subsequently, you can observe the harmonics of the string one by one by pressing ▲ *Up*.

(After each frequency jump: Wait until the previous resonance has died out and the new one has evolved.)

## Sweep

The function generator can perform linear and logarithmic frequency sweeps.

(A linear sweep passes "equal numbers of Hz per second", a logarithmic one "equal numbers of octaves per second").

A sweep is defined by four parameters: *type* (Lin/Log), *start frequency*, *stop frequency*, and *sweep time*. To set up these parameters in the order mentioned, you press *Sweep* and use the knob *Frequency*. The display indicates the parameter currently being set.

The step function can be used when setting start and stop frequencies.

After the four parameters have been set, the next press of *Sweep* activates the sweep.

The sweep function is interrupted by yet another press of *Sweep*.

While a sweep is active, the frequency span may be changed with the buttons ▲ *Up* and ▼ *Down*. The start frequency changes as described in the section *The step function*. The stop frequency is changed according to the following principle: For a linear sweep the difference *stop frequency* – *start frequency* is preserved. For a logarithmic sweep the fraction *stop frequency* / *start frequency* is preserved.

## Application example – Sweep

To demonstrate the sharp resonances in Chladni plates, sweep the frequency slowly through an interval. With a little sand on the plate, the nodal patterns will stand out.

Press *Sweep* and rotate the (frequency-) knob slightly until the display shows *Lin*.

Press again (*Start* is now blinking in the display), and set the desired start frequency. (50 Hz is a good bet.)

Press *Sweep* again (*Stop* blinking) and set the desired stop frequency. (600 Hz is a good bet.)

After the next push (*Time* flashes on the display) sweep time is set to for example 20 seconds.

Now the sweep is ready and will start running at the next press of *Sweep*.

Sweep parameters are remembered. If a single parameter is to be changed, simply step forward by pressing the *Sweep* button. The values not actively changed remain the same.

## Advanced use

The function generator behaves as a 1 MB memory stick when connected to a PC via USB.

Note that it is very important that you close the USB connection properly down before the cable is removed, or any of the appliances turned off. Always use the function "Safely remove hardware" when you are finished using the connection.

The drive contains the following folders:

- The root folder "**"**"  
(may eventually contain an editable setup file)
- "firmware"**
- " waveforms\predefined"**  
(hidden folders – avoid changing anything here)
- "settings"** (hidden folder – editable text files )
- "waveforms\Custom"** (the definition of a custom-defined waveform may be placed here)

## Automatic mode

It is possible to define a sequence of settings to be carried out by the function generator. This may just be the automation of a single setting - or it could be a long series of e.g. frequency changes.

Setting up automatic mode is done in the file **"setup.zup"**. The file should be in the USB drive's root folder. The possible commands are specified below.

Activation of the automatic mode is achieved by holding the Display button down and pressing the *f<sub>0</sub>/Mode*. The function is indicated as AM in the display.

The generator returns to normal operation by pressing a button or by turning the frequency knob.

It is recommended that you take a backup of the drive before making changes.

## Commands in "setup.zup"

This text file contains one command per line. Each command line starts with a letter, eventually followed by a number. Empty lines and comments are allowed. A comment is a line that does not start with any of the letters below. It is recommended always to start a comment with two forward slashes: *"//"*.

The commands are executed from top to bottom (except when an R-command is executed - see below).

D	Dump	Prints the current settings to the file dump.txt. Each line consists of three fields: Parameter name e.g. "Waveform" Parameter value e.g. "Ramp Up." Corresponding command e.g. "W6"
Fx	Freq.	Set the frequency to the value x. For example F234500 will set f to 234.5 kHz.
Wn	Waveform	The number n indicates the waveform. W1 means sine, W6 is upward ramp. If there is no custom waveform, W8 means the same as W1.
Bx	Base freq.	Sets $f_0$ to the frequency x.
Mn	Step Mode	M1: harmonics, M2: octaves, M3 decades.
Ni	N-value	The current N-value is set to the integer i – which can take both positive and negative values. If step mode is M1, i=0 and i=-1 will be ignored
Sn	Sweep mode	S0: normal operation, S1: linear sweep, S3: logarithmic sweep.
Ix	Initial freq.	Set the start frequency of sweep to x.
Ex	End freq.	Set the stop frequency of the sweep to x.
Tx	Sweep Time	The time x is a decimal number in seconds.
Px	Pause	Wait x seconds before the next line is executed. The signal from the generator continues with the current parameters.
R	Repeat	Jump to the top of the file and repeat from there. Lines following an R-command are never executed.

Frequencies can be specified with 0.001 Hz resolution. Precision is better than 0.005 %.

## Custom-defined waveform

There are two possible formats for the specification of a waveform. The file is placed in the folder "**\waveforms\Custom**".

If you want the display to show the amplitude (peak value) correctly, you must ensure that the maximum value occurs in the file (4095 for "**custom.csv**" resp. 100 for "**custom.wdf**").

### "custom.csv" (comma separated values)

This text file consists of 16384 lines, each consisting of one integer between 0 and 4095 (incl.).

The value 0 corresponds to the voltage -100 % and the value 4095 corresponds to 100 %.

Comments may be inserted in the form of lines that do not start with a digit.

The content is generated easily by using a spreadsheet – copy a column to a text file and rename.

If "**custom.csv**" exists, the function generator creates the file "**custom.wfr**" in its internal format. It is a binary format and is not directly editable. This transformation takes time, so the first time you select the waveform *Prog.* you will experience a few seconds response delay. Then "**custom.csv**" is automatically renamed "**custom. \$cs**".

### "custom.wdf" (waveform description file)

This file is a quick way to define a waveform that consists of line segments. Note that if "**custom.wfr**" exists, "**custom.wdf**" is ignored.

Each line in the file describes an endpoint of a linear ramp given by two numbers:

A **time** (0 to 100 %) and a **voltage** (from -100 % to 100 %) - separated by semicolons or tabulation.

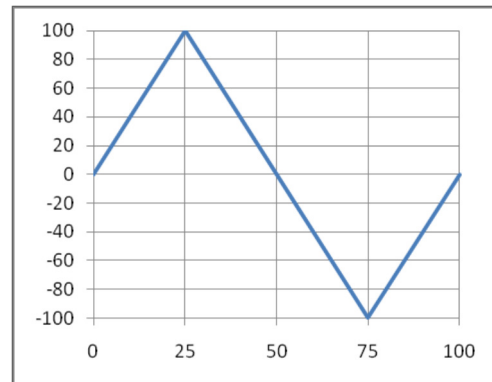
If the time entry in the first point is not 0, the curve starts at (0,0). If the last time is less than 100, the last defined voltage continues for the rest of the period.

Comments may be inserted in the form of lines that do not start with a digit

### Waveform description file – example

```
//Triangle Wave
25; 100
75; -100
100; 0
```

(Since the first time is 25 %, the starting voltage is implicitly 0 V.)

Note that it is very important that you close the USB connection properly down **before** the cable is removed, or any of the appliances turned off.

**Always** use the function "Safely remove hardware" when you are finished using the connection.

## System parameters

The file "**system.ini**" lies in the hidden folder "**settings**". This file makes it possible to specify a few more unusual parameters. Here we shall only mention a few. Normally there will be no need to change these parameters:

*VoltComp = ON/OFF*

Analogue components of the generator have a finite bandwidth, and the amplitude therefore decrease a bit at high frequencies. The decrease depends on the waveform. With VoltComp = ON, the amplitude, as shown in the display, reflect this. With VoltComp = OFF the "ideal" value is displayed.

*Encoder = 0/1/2/3/4*

The frequency control is as previously mentioned speed sensitive. You can choose from 5 different built-in acceleration characteristics.

(This parameter will be ignored if there is a file called "**encoder.ini**" in the folder.)

Note - In the "**system.ini**" file there is one space on each side of the equal sign.

## Specifications

### Waveforms

Bipolar:	Sine, triangle, square
Positive:	Square pulse, triangle pulse, ramp up, ramp down
Additionally user defined, arbitrary waveform.	
Distortion (sine):	< 0.1 % up to 20 kHz; < 1 % otherwise

### Frequency range

50 $\Omega$ output:	0.001 Hz to 10.00 MHz
Sync output:	0.001 Hz to 10.00 MHz
Power output:	0.001 Hz to 100.0 kHz
Frequency stability:	Better than 0.005 %

### Amplitude

50 $\Omega$ output, no load:	0 to 10 V (20 V p-p for bipolar waveforms)
50 $\Omega$ output, 50 $\Omega$ termination:	0 to 5 V (10 V p-p for bipolar waveforms)
Sync output:	5 V (TTL signal: 0 to 5 V)
Power output:	0 to 10 V (20 V p-p for bipolar waveforms)

### Max current

50 $\Omega$ output, short circuit:	200 mA (only briefly)
50 $\Omega$ output, 50 $\Omega$ termination:	100 mA (unlimited)
Sync output, short circuit:	100 mA (unlimited)
Power output:	1 A

### Misc.

Power consumption:	85 W (max) ; 21 W (idle)
Dimensions (W x D x H):	312 x 205 x 117 mm

### Mains voltage

	Model 250350	Model 250351
Mains voltage:	230 V ; 50-60 Hz	115 V ; 50-60 Hz
Fuse:	1 A (slow)	2 A (slow)

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© Frederiksen Scientific A/S