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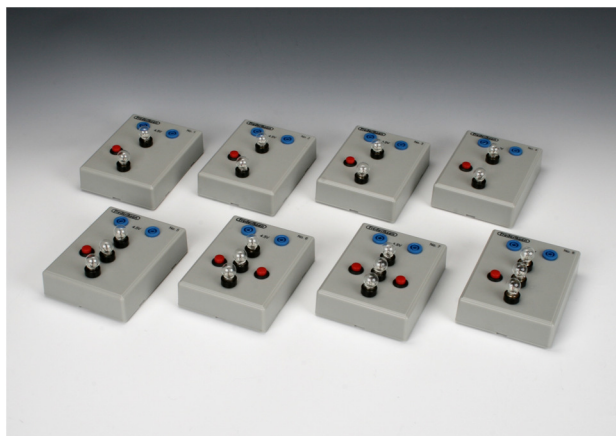
Eight different hidden circuits with switches and incandescent lamps.

Each box is powered by a 4.5 V battery or a power supply (not included).

The circuits can be investigated by pushing the switch buttons and by screwing the bulbs out or in.

The students are challenged to determine the configuration of the circuits and to draw the schematics diagrams for each of them.

There is some progression from box No. 1 to No. 8.

Relevant documents

A corresponding **lab manual** for the students can be downloaded from www.frederiksen.eu – search for item number 429800 or the manual number **136040**).

A **teacher's guide** is included which among other things contains solutions (**not** available for download).

Equipment required

The circuits can be powered from a 4.5 V battery:

350555 Battery, 3R12 4,5 V

Use lab leads and a pair of crocodile clips to make contact with the battery.

As an alternative, a power supply can be used. Almost any unit will work: for instance 361600, 261700, 361055, 361870, 364000 – as long as it can deliver 4 V and 0.5 A AC or DC. When using a power supply, it is best not to go all the way up to 4.5 V.

Spare parts

The light bulbs are of the following type:

425015 Bulb, 3.5 V, 0.2 A (Qty. 10)

A sufficient number of bulbs are included but it is recommended to have some spare ones ready.

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© Frederiksen Scientific A/S