


LabField

067960 Roller Mixer, 6 rollers

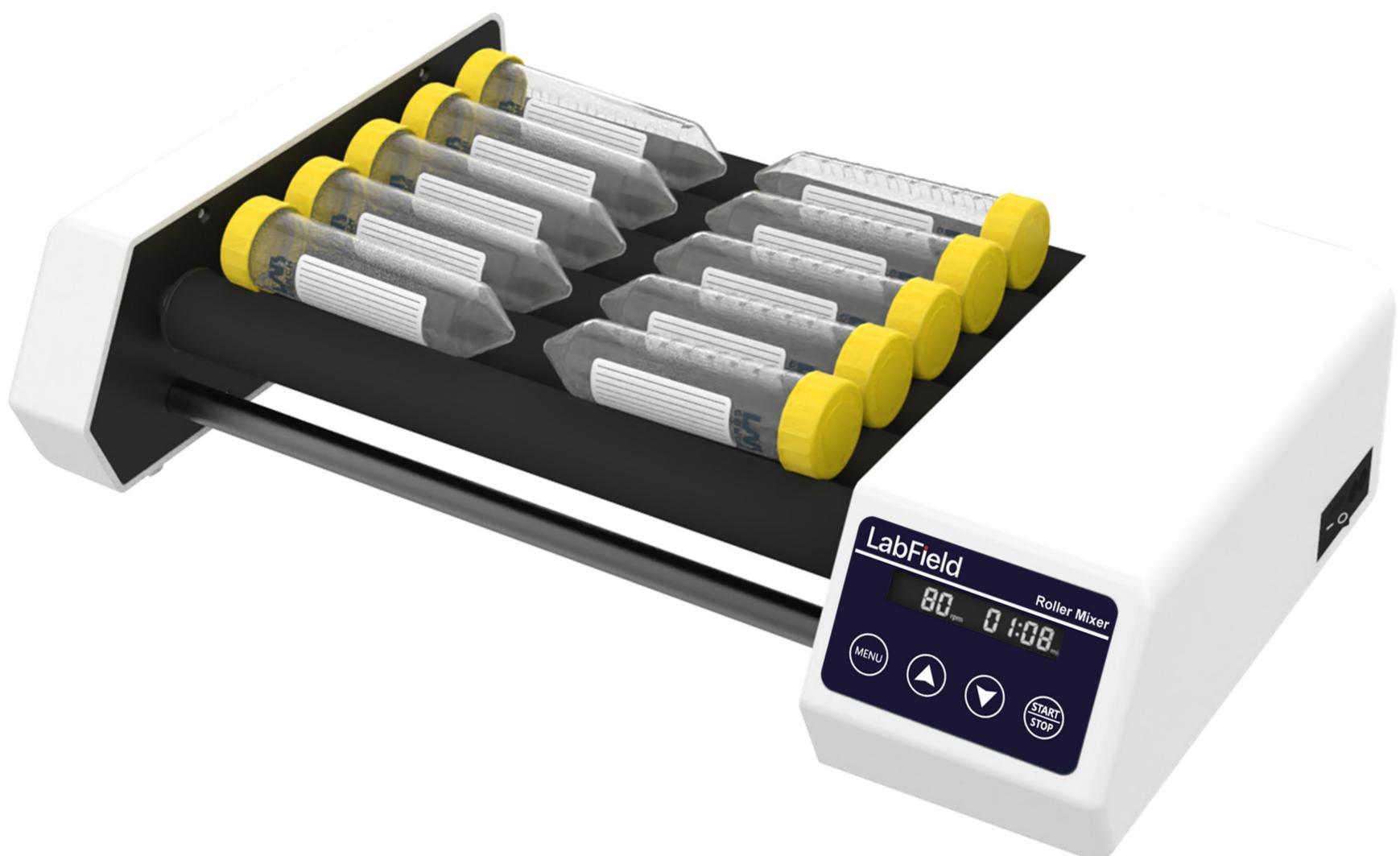


User manual


I: Product brief

Labfield mixers are designed and manufactured according to international advanced design concepts and technologies. The mixers are highly efficient and gentle and are suitable for different experiments such as blood analysis and sample mixing.

II: Product overview

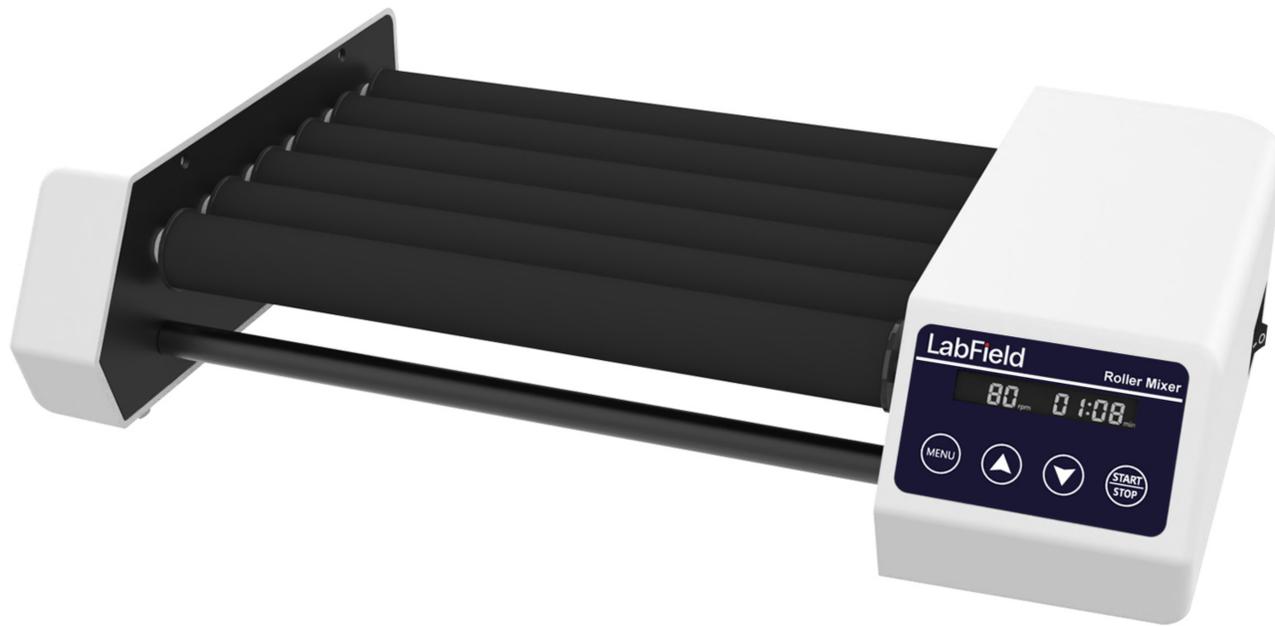

III: Product introduction

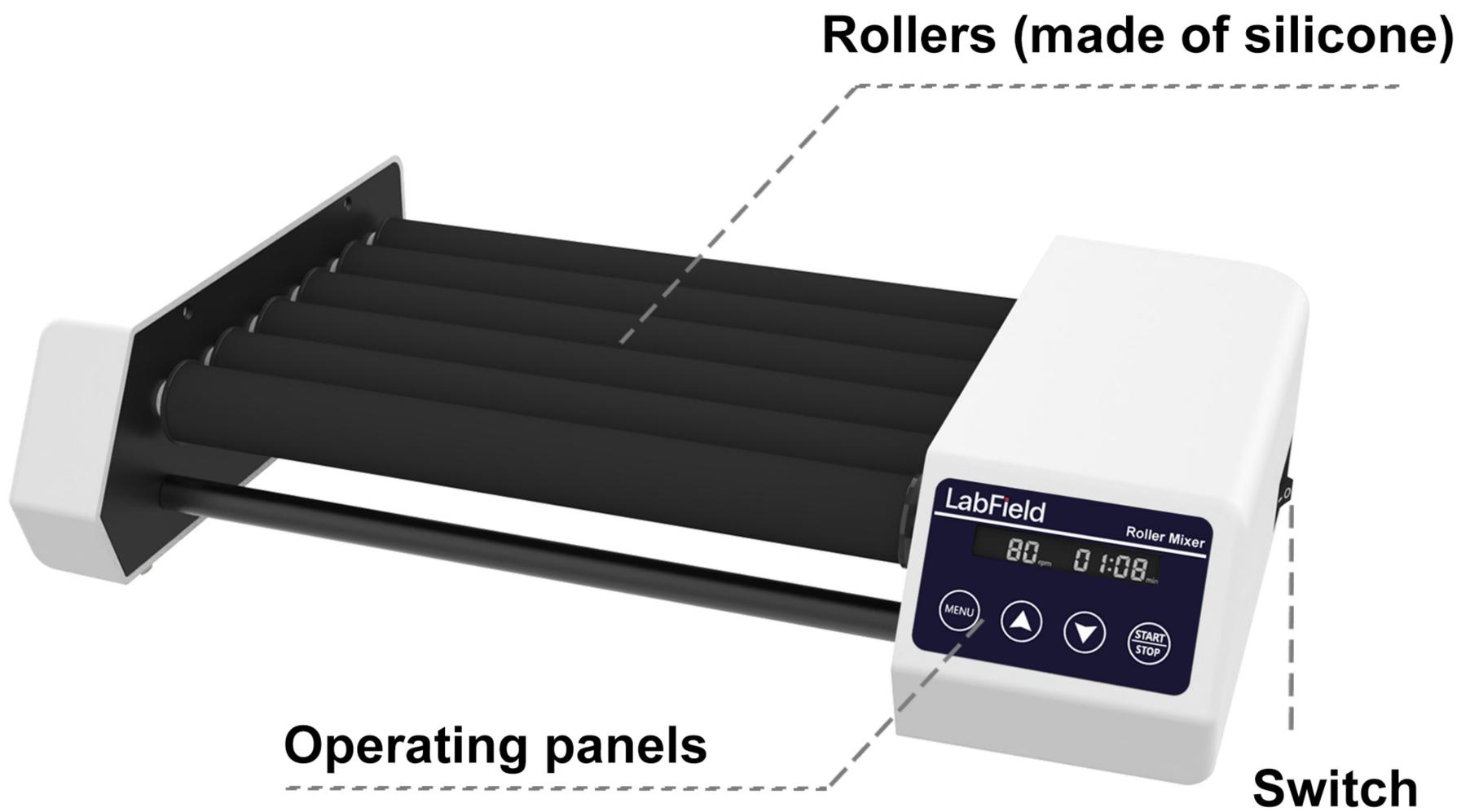

IV: Product data

Type	Horizontal rolling
Rollers	6 rollers
Max Capacity	10*50ml Centrifuge tube / 0.5kg
Control mode	Digital Key
Speed	20-80 rpm
Timing	1min - 23h59min
SPECS	φ6-34mm
Power	25W
Voltage	AC110-220V
Dimension(mm)	440*255*105

V: Correct usage

Application

- Suitable for mixing operations in different applications: electron microscopy analysis, blood cell suspensions, homogeneous solid blood and plasma, tissue specimens, washing precipitates, disintegration sheet analysis, extraction process, dissolution slow reaction, rotary sealing, etc.

Usage

- Widely used in scientific research institutions, medical schools, health care and other units, molecular biology, virology, microbiology, pathology, immunology and other laboratories.

Steps for usage

- Plug in the power supply and turn on the switch;
- Place the sample tube/reagent bottle on the roller;
- MENU key to switch time/speed, up and down keys to adjust parameters;
- Press the START/STOP keys to start/stop the device.

Instrument using environment

- Indoor use
- Altitude: $\leq 2000\text{m}$
- The instrument is suitable for ambient temperature range of $+5^{\circ}\text{C}\sim+40^{\circ}\text{C}$
- Instrument operation suitable for relative humidity range of $\leq 80\%$
- No vibrations and airflow affecting performance are present
- There are no conductive dust, explosive gases and corrosive gases in the surrounding air
- Keep a safety distance of 10cm between the front and back left and right when the instrument is working

Safety Precautions

- It is strictly forbidden to unplug the power connector and toggle the switching power button when your hands are stained with liquid!
- It is strictly forbidden to unplug the power plug when the instrument is live!
- It is strictly forbidden to maintain and scrub the instrument when it is powered on!
- It is strictly forbidden to install the instrument on an uneven, shaky and shaky work surface!

Delivery list

- Host
- Power cable
- User manual

